



Product Specifications

Electrophoresis Reagents, Polymerase Chain Reaction
Custom Primers and Probes
Hybridization and Detection Reagents

PCR Additives

Store at -20°C

	Catalog Number	Description	Unit Size
<input type="checkbox"/>	40-3020-05	dNTP 10mM	0.5 ml
<input type="checkbox"/>	40-3021-11	dNTP 2mM (10X)	1.1 ml
<input type="checkbox"/>	40-3022-16	MgCl ₂ ; 25 mM	1.6 ml
<input type="checkbox"/>	40-3023-16	MgCl ₂ ; 50 mM	1.6 ml
<input type="checkbox"/>	40-3031-10	DMSO	1 ml
<input type="checkbox"/>	40-3053-10	TMAC (Tetramethyl ammonium chloride) 100 mM	1 ml
<input type="checkbox"/>	40-3059-10	KCl 300 mM	1 ml
<input type="checkbox"/>	40-3032-10	Betaine; 5M	1 ml

PCR Additives

DNA polymerases need to elongate rapidly and accurately to function effectively *in vivo* and *in vitro*, yet certain DNA regions appear to interfere with their progress. One common problem is pause sites, at which DNA polymerase molecules cease elongation for varying lengths of time. Many strong DNA polymerase pauses are at the beginnings of regions of strong secondary structure such as template hairpins (1). Taq polymerase use in PCR suffers the same fate and GC-rich DNA sequences often require laborious work to optimize the amplification assay. The GC-rich sequences possess high thermal and structural stability, presumably because the high duplex melting temperature that permits stable secondary structures to form, thus preventing completion of a faithful replication (2).

Nucleotide analog 7-deaza dGTP is effective in reducing the secondary structure associated with GC rich region by reducing the duplex stability(4). Betaine, DMSO and formamide reduces the T_m and the complex secondary structure thus the duplex stability (1-5). Tetramethyl ammonium chloride (TMAC) actually increases the specificity of hybridization and increases the T_m. The use of TMAC is recommended in PCR conditions using degenerate primers.

These PCR additives and enhancing agents have been used to increase the yield, specificity and consistency of PCR reactions. These additives may have beneficial effects on some amplification and it is impossible to predict which agents will be useful in a particular context and therefore they must be empirically tested for each combination of template and primers.

PCR Additives		
Additive	Purpose & Function	Concentration
7-deaza-2'-deoxyguanosine; 7-deaza dGTP	GC rich region amplification. Reduce the stability of duplex DNA	Totally replace dGTP with 7-deaza dGTP; or use 7-deaza dGTP: dGTP at 3:1
Betaine (N,N,N-trimethylglycine = [carboxymethyl]trimethylammonium)	Reduces T _m facilitating GC rich region amplification. Reduces duplex stability	Use 3.5M to 0.1M betaine. Be sure to use Betaine or Betaine (mono)hydrate and not Betaine HCl.
BSA (bovine serum albumin)	BSA has proven particularly useful when attempting to amplify ancient DNA or templates which contain PCR inhibitors such as melanin.	BSA concentration of 0.01µg/µl to 0.1µg/ µl can be used.
DMSO (dimethyl sulfoxide)	DMSO is thought to reduce secondary structure and is particularly useful for GC rich templates.	DMSO at 2-10% may be necessary for amplification of some templates, however 10% DMSO can reduce <i>Taq</i> polymerase activity by up to 50% so it should not be used routinely.
Formamide	Reduces secondary structure and is particularly useful for GC rich templates.	Formamide is generally used at 1-5%. Do not exceed 10%.
Non-ionic detergents e.g. Triton X-100, Tween 20 or Nonidet P-40 (NP-40)	Non-ionic detergents stabilise <i>Taq</i> polymerase and may also suppress the formation of secondary structure.	0.1-1% Triton X-100, Tween 20 or NP-40 may increase yield but may also increase non-specific amplification. As little as 0.01% SDS contamination of the template DNA (left-over from the extraction procedure) can inhibit PCR by reducing <i>Taq</i> polymerase activity to as low as 10%, however, inclusion of 0.5% Tween-20 or -40 will effectively neutralize this effect.
TMAC (tetramethylammonium chloride)	TMAC is used to reduce potential DNA-RNA mismatch and improve the stringency of hybridization reactions. It increases T _m and minimizes mis-pairing.	TMAC is generally used at a final concentration of 15-100mM to eliminate non-specific priming.

Appendix

PCR Components and Analysis

PCR buffer conditions vary and it is imperative to optimize buffer conditions for each amplification reaction. At Gene Link most amplification reactions have been optimized to work with the following standard buffer condition, unless indicated.

Standard Gene Link PCR Buffer	
10 X PCR buffer	1 X PCR buffer
100 mM Tris-HCl pH 8.3	10 mM
500 mM KCl	50 mM
15 mM MgCl ₂	1.5 mM
0.01% Gelatin	0.001%

dNTP Concentration

Standard dNTP concentration of 0.2mM of each base is used. See section on PCR additives when dNTP concentration is changed.

Recipe	
2.0 mM dNTP Stock Solution Preparation*	
100 mM dGTP	100 µl
100 mM dATP	100 µl
100 mM dTTP	100 µl
100 mM dCTP	100 µl
Water	4.6 ml
Total Volume	5 ml
*Aliquot and freeze	

MgCl₂ Concentration

The concentration of Mg⁺⁺ will vary from 1-5 mM, depending upon primers and substrate. Since Mg²⁺ ions form complexes with dNTPs, primers and DNA templates, the optimal concentration of MgCl₂ has to be selected for each experiment. Low Mg²⁺ ions result in a low yield of PCR product, and high concentrations increase the yield of non-specific products and promote mis-incorporation. Lower Mg²⁺ concentrations are desirable when fidelity of DNA synthesis is critical. The recommended range of MgCl₂ concentration is 1-4 mM, under the standard reaction conditions specified. At Gene Link, using the standard PCR buffer with KCl, a final dNTP concentration of 0.2 mM, a MgCl₂ concentration of 1.5 is used in most cases. If the DNA samples contain EDTA or other chelators, the MgCl₂ concentration in the reaction mixture should be raised proportionally. Given below is an MgCl₂ concentration calculation and addition table using a stock solution of 25 mM MgCl₂.

MgCl ₂ Concentration & Addition Table								
Final concentration of MgCl ₂ in 50µl reaction mix, (mM)	1.0	1.25	1.5	1.75	2.0	2.5	3.0	4.0
Volume of 25mM MgCl ₂ , µl	2	2.5	3	3.5	4	5	6	8

Primer Concentration

The final concentration of primers in a PCR reaction is usually 0.5 to 1 μM (micromolar). This is equivalent to 0.5 to 1 pmol/ μl . For a 100 μl reaction you would add 50 to 100 pmols. At Gene Link we use 0.5 pmol/ μl in the final PCR.

Primer Reconstitution

Stock Primer Mix: Dissolve the supplied 10 nmols of lyophilized Genemer™ in 100 μl sterile TE. The 10 nmols of primers when dissolved in 100 μl will give a solution of 100 μM i.e. 100 pmols/ μl .

Primer Mix: Prepare a 10 pmols/ μl Primer Mix solution by a ten fold dilution of the stock primer mix.

Example: Add 180 μl sterile TE to a new tube, to this tube add 20 μl of primer stock solution. Label this tube as Primer Mix 10 pmols/ μl .

Amplification Thermal Cycling

Hot Start: It is essential to have a 'Hot Start' profile for amplification of any fragment from a complex template like human genomic DNA. Taq polymerase has low activity at room temperature and it is essential to minimize any mis-priming in the first cycle of amplification. A typical hot start profile is given below. Various enzyme preparations are available which are activated by heat in the first cycle. A simple hot start protocol is given below that can be used with regular Taq polymerase. See the section on PCR additives for amplification of products from high GC content templates.

Hot Start		
Step	Time & Temperature	Cycles
Initial Denaturation	95 °C for 5 minutes	1
Annealing	60 °C Hold Infinity	Hold
Comments: Add Taq premix while on hold.		

Amplification File

The initial denaturation step at 94 °C for 30 seconds is sufficient for all templates. The number of cycles is usually set to 30 and is sufficient to amplify 1-10 μg of product depending on the initial concentration of template. A higher number of cycles from 35-45 cycles may be used, but internal priming on the product and over amplification of unwanted bands often result from over-cycling. Generally, it is better to focus on optimizing reaction conditions than to go beyond 35 cycles.

Typical Amplification File			
Step	Temperature	Time	Cycles
Denaturation	94 °C	30 sec.	30
Annealing	*	30 sec.	
Elongation	72 °C	30 sec.	
Fill in Extension	72 °C	7 minutes	1
Hold	4 °C	Infinity	Hold
*Based on the Tm of the primers. Usually varies from 50°C to 65°C			

Typical Reaction Premix

Typical PCR Premix (/50µl)	
Component	Volume
10 x PCR Buffer	5 µl
2.0 mM dNTP mix (each)	5 µl
Primer Mix (10 pmol/µl each) or 2.5µl of 10 pmol/µl of individual primer (final 25 pmol of each primer/50µl)	2.5 µl
H ₂ O	37.5 µl
Total Volume	50 µl

Typical PCR Reaction Mix

PCR reaction (/50µl)	
Component	Volume
PCR premix	45 µl
100ng/µl diluted DNA	1 µl
Hot start and then add	
Taq premix	5 µl


Taq Premix Preparation

Taq Premix (/50µl)	
Component	Volume
PCR Premix	6 µl
Taq polymerase (5 u/µl)	0.25µl
Add 5µl/50 µl rxn. After initial denaturation	
Use 2.5 units of Taq for 100 µl reactions. Taq is usually supplied at a concentration of 5 units/µl	

Yield and Kinetics

The target will be amplified by up to 10⁶ fold in a successful reaction, but the amplification will usually plateau at 1-10µg. Thus, 1 pg of target sequence in the reaction is a good place to begin.

PCR reactions produce product in a nonlinear pattern. Amplification follows a typical exponential curve until some saturation point is reached. Generally products will not be further amplified once 1-5 µg has been generated. Saturation by one product of a reaction does not always prevent further amplification of other generally unwanted products. Over-cycling may decrease the quality of an otherwise good reaction. When first optimizing a reaction, it is advisable to take samples every 5 or 10 cycles to determine the number of cycles actually needed.

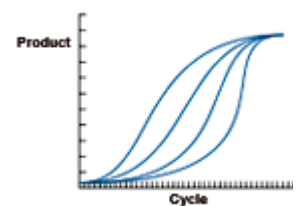

References:

1. Kovarova, M; and Draber, P; (2000) New Specificity and yield enhancer for polymerase chain reactions (2000) Nucl. Acids. Res. 28: e70.
2. Henke, W., Herdel, K., Jung, K., Schnorr, D. and Stefan A. Loening, S. (1997) Betaine improves the PCR amplification of GC-rich DNA sequences. Nucl. Acids Res. 25: 3957-3958.
3. Daniel S. Mytelka, D.S., and Chamberlin, M.J.,(1996) Analysis and suppression of DNA polymerasepauses associated with a trinucleotide consensus. Nuc. Acids Res.,. 24:2774–278.
4. Keith, J. M., Cochran, D.A.E., Lala, G.H., Adams, P., Bryant, D.and Mitchelson, K.R. (2004) Unlocking hidden genomic sequence. Nucl. Acids Res. 32: e35.
5. Owczarzy, R., Dunietz, I., Behlke, M.A., Klotz, I.M. and Joseph A. Walder. (2003) Thermodynamic treatment of oligonucleotide duplex–simplex equilibria. PNAS, 100:14840-14845.

Ordering Information

Product	Catalog No.	Unit Size
Taq DNA Polymerase; 400 units; 5 µl/µl; 80 µl	40-5200-40	400 units
Taq PCR Kit; 200 x 50 µl reactions	40-5211-01	200 reactions
Taq PCR Kit with controls; 200 reactions	40-5212-01	200 reactions
PCR Master Mix (2X); 100 x 50 µl reactions (2 tubes x 1.3 ml)	40-5213-01	100 reactions
PCR Master Mix (2X); 200 x 50 µl reactions (4 tubes x 1.3 ml)	40-5213-02	200 reactions

Related Products Ordering Information

PCR Reagents

Product	Catalog No.	Unit Size
Taq DNA Polymerase 300 units; 5 µl/µl; 60 µl	40-5200-30	300 units
PCR Buffer Standard (10 X)	40-3060-16	1.6 ml
PCR Buffer Mg Free (10 X)	40-3061-16	1.6 ml
Taq Polymerase Dilution Buffer; 1 ml	40-3070-10	1 ml
dNTP 2mM (10X)	40-3021-11	1.1 ml
MgCl ₂ ; 25 mM	40-3022-16	1.6 ml
Omni-Marker™ Universal Unlabeled	40-3005-01	100 µl
Primer and Template Mix; 500 bp; 40 reactions	40-2026-60PT	100 µl
Nuclease Free Water	40-3001-16	1.6 ml
DMSO	40-3031-10	1 ml
TMAC (Tetramethyl ammonium chloride) 100 mM	40-3053-10	1 ml
KCl 300 mM	40-3059-10	1 ml
Betaine; 5M	40-3032-10	1 ml

Omni-Marker™

Product	Catalog No.	Unit Size*
Omni-Marker™ Universal unlabeled	40-3005-01	100 µl
Omni- Marker™ Universal unlabeled	40-3005-05	500 µl
Omni-Marker™ Universal unlabeled	40-3005-10	1 ml
Omni- Marker™ Low unlabeled	40-3006-01	100 µl
Omni-Marker™ Low unlabeled	40-3006-05	500 µl
Omni- Marker™ Low unlabeled	40-3006-10	1 ml
Omni-Marker™ GScan-2 Tamra labeled 50 bp - 600 bp	40-3062-01	100 µl
Omni-Marker™ GScan-2 Tamra labeled 50 bp - 600 bp	40-3062-05	500 µl

Buffers & Reagents

Product	Catalog No.	Unit Size
Agarose Tablets, 0.5 gm each	40-3011-10	100 tablets
Agarose LE Molecular Biology Grade; 100 gms	40-3010-10	100 g
Agarose LE Molecular Biology Grade; 500 gms	40-3010-50	500 g
Hybwash A, Hybridization Wash Solution	40-5020-20	200 ml
Hybwash B, Hybridization Wash Solution	40-5021-10	100 ml
TAE Buffer; 50X Concentrate; 100 ml	40-3007-01	100 ml
TAE Buffer; 50X Concentrate; 1000 ml	40-3007-10	1000 ml
TBE Buffer; 5X Concentrate	40-3008-10	1000 ml
10x Washing buffer	40-5025-20	200 ml
10% Blocking solution	40-5026-10	100 ml
Seq. Loading buffer	40-5027-00	1 ml
10x AP Detection buffer	40-5031-10	100 ml
Lumisol™ I Hybridization Solution; contains formamide	40-5022-20	200 ml
Lumisol™ II Hybridization Solution; for non-toxic hybridizations	40-5023-20	200 ml
Lumisol™ III Hybridization Solution; for oligo probes	40-5024-20	200 ml

Loading Buffers

Product	Catalog No.	Unit Size
Loading Buffer 5X BPB/XC non-denaturing	40-3002-01	100 µl
Loading Buffer 5X BPB/XC non-denaturing	40-3002-10	1 ml
Loading Buffer 5X Orange G/XC non-denaturing	40-3004-01	100 µl
Loading Buffer 5X Orange G/XC non-denaturing	40-3004-10	1 ml
Loading Buffer 2X BPB/XC Denaturing for Sequencing	40-5027-01	100 µl
Loading Buffer 2X BPB/XC Denaturing for Sequencing	40-5027-10	1 ml

Prices subject to change without notice

All Gene Link products are for research use only

Gene Link Distributors

Argentina

Migliore Laclaustra S.R.L.
Buenos Aires 1055 Argentina
(11) 4372 9045
(11) 4373 8052
Email :
miglac@migliorelaclaustra.com.ar

Belgium

Gentaur
Kampenhout 1910 Belgium
0032 16 58 90 45
0032 16 50 90 45
Email : info@gentaur.com
Web : www.gentaur.com

Brazil

LGC do Brasil
Rio de Janeiro 20770-270 Brazil
Tel 21-592-6642
Toll Free 0800- 407047
Fax 21-2593-3232
Web : www.lgcscientific.com

China

Hangzhou AOWEI Bioengineering
Hangzhou China
Tel: 86-871-56773688 /
86-571-56773688
86-571-56773066
Email : liuwg@aowei21.com
Web : www.aowei21.com

Egypt

Lab Technology
P.O.Box 5959 Heliopolis West
Cairo 11351 Egypt
20-2-2622 6761
20-2-2622 6760
Email : subhi@labtechnology.net

India

Lab Equipments & Chemicals
Kolkata 700 001 India
Tel: 91-33-2235-1958/1806/0274
Fax: 91-33-2221-5682
Email : info@labquipindia.com
Web : www.labquipindia.com

Ireland

Isis Ltd.
Bray, Co Wicklow Ireland
Tel: 00353-1-2867777
Fax: 00353-1-2867766
Email : enquiries@isisco.ie
Web :
www.isisco.ie/isis/Main/Home.htm

Italy

DBA Italia S.R.L
Milano Italy
Tel: 39-02-26922300
Fax: 39-02-26923535
Email : info@dbaitalia.it
Web : www.dbaitalia.it

Japan

Funakoshi Co., Ltd.
Tokyo 113-0033 Japan
Tel: (3) 5684-1620
Fax: (3) 5684-1775
Email : reagent@funakoshi.co.jp
Web : www.funakoshi.co.jp

Jordan

LeanGene for Medical & Electronic
Supplies
Amman 11110 Jordan
(962) 6 46 55 401
(962) 6 46 55 402
Email : mgr@leangene-jo.com

Jordan

Advanced Scientific Consulting
Group Co.
City Terminal bldg. Office 303
[144753]
Amman Jordan
962 6 5812210
962 6 5812214
Email : asem@advanced.com.jo
Web : www.advanced.com.jo

Jordan

CGenomix
Amman Jordan
+ 962 (6) 55 22 365
+ 962 (6) 55 22 367
Email : info@cgenomix.com
Web : www.cgenomix.com

Korea

Genes Laboratories Inc.
Soojeong-gu, Seongnam-si
(461-701)
Gyeonggi-do Korea
82-(0)31-889-3085
82-(0)31-696-5453
Email : info@geneslabs.com
Web : www.geneslabs.com

Korea

GenomicBase Co. Ltd.
Dongdaemun-Gu [130-842]
Seoul Korea
Tel: 82. 502.335.4926
Fax: 82. 502.335.4927
Email : info@genomicbase.com
Web : www.genomicbase.com

Kuwait

Tri Alpha Co.
Al-Awadi Tower #3, 6th Floor
Kuwait City 72263 Kuwait
(00965) 22 43 13 44
(00965) 22 43 51 06
Email : walter@trialphaco.com

Latvia

Valdai
Latvia LV-1010 Latvia
Tel: + 371-7227620
Fax: + 371-7326091
Email : Valdai@apollo.lv

Lebanon

DPC/Lebanon SARL
PO BOX 55440
Beirut Lebanon
Tel: (961) 502 812
Fax: (961) 805 851
Email : dpcliban@dm.net.lb
Web : www.dpclub.com

Malaysia

Nano Life Quest Sdn Bhd
47100 Puchong, Selangor D. E.
MALAYSIA
Tel : +603 8062 4678
Fax : +603 8062 4678
Email : info@nanolifequest.com
Web : www.nanolifequest.com

Norway

AME Bioscience A/S
Torsoed 3135 Norway
47-3361-0718
47-3340-1444
Email: aewing@amebioscience.com

Pakistan

Worldwide Scientific
30-Ferozepur Road
Lahore 54000 Punjab Pakistan
Tel: 92-42-755-2355
Fax: 92-42-755-3255
Email : wws@brain.net.pk

Saudi Arabia

Labs Care
Riyadh 11322 Saudi Arabia
Tel: (1) 465-9452
Fax: (1) 465-9452
Email : manager@labscare.com
Web : www.labscare.com

Singapore

Sciencewerke Pte Ltd
#04-21 S139550
Singapore
Tel: +65 777 1045
Fax: +65 777 3054
Email : enquiry@sciencewerke.com
Web : www.sciencewerke.com

Spain

bioNova científica, s.l.
28007
Madrid Spain
Tel: +34 91 551 54 03
Fax: +34 91 433 45 45
Email : info@bionova.es
Web : www.bionova.es

Spain

Attendbio Research SL
08018
Barcelona Spain
Tel: +34 934019842
Fax: +34 934019843
Email : attendbio@attendbio.com
Web : www.attendbio.com

Taiwan

Watson Biotechnology Co., Ltd.
Xinzhuang Dist.
New Taipei City 242 Taiwan
Tel: 886-2-89916881
Fax: 886-2-89936967
Email : tensci.gene@msa.hinet.net
Web : www.tenscigene.com.tw

Thailand

Qualitech Instruments Co.,Ltd
Ladprow
Bangkok 10230 Thailand
Tel: (662) 538-1843
Fax: (21) 2593-3232
Email : sales@qualitech-instru-
ments.com
Web : www.qualitech-
instruments.com

Turkey

Diagen Biyoteknolojik Sistemler
Halk Sokak 27/2 Sihhye
Ankara 06420 Turkey
Tel: (312) 432-0611
Fax: (312) 432-3595
Email : info@diagen.com.tr
Web : www.diagen.com.tr

UK & Switzerland

Molecular Solutions Europe Ltd.
London EC1V 4PY UK UK &
Switzerland
Tel: +44-(0)870-1995067
Fax: +44-(0)207-2539040
Email : info@mseu.co.uk
Web : www.mseu.co.uk

United Arab Emirates

Al Nawras Medi-Lab Supplies
Industrial Area No. 11
Sharjah United Arab Emirates
Tel: (9716) 5344550
Email : alnawras@eim.ae

Gene Link, Inc.

190 Saw Mill River Road
Hawthorne, NY 10532
USA

Telephone:

1-800-GENE-LINK
(914) 769-1192

Facsimile:

1-888-GENE-LINK
(914) 769-1193

Email:

CUSTOMER SERVICE

cust_service@genelink.com

SALES

sales@genelink.com

CUSTOM OLIGO ORDERS

oligos@genelink.com

TECHNICAL SUPPORT

support@genelink.com

SEQUENCING/GENOTYPING

genotyping@genelink.com

ALL ORDERS

orders@genelink.com



Gene Link™