



Certificate of Analysis & Product Manual

SmartBase™ siRNA Custom siRNA duplex,
siRNA fluorescent and modified
Guaranteed RNAi Explorer

siRNA & RNAi Explorer™

Catalog No.: See Material Supplied List

Storage Condition: -20°C

For Research Use Only. Not for use in diagnostic procedures for clinical purposes



Certificate of Analysis & Product Specifications

Please consult and save enclosed 'Order Specifications' sheet with a clearly marked order number and or lot number, this gives order specific data including the exact nmol quantity supplied of each duplex.


Product Label Information

Guaranteed RNAi Explorer™

	Content	Catalog No.	Product	Size
REF	<input type="checkbox"/>	27-6402-02	3 Target siRNA's & 1 negative control duplex. Standard purity, 2 nmol each	2 nmols
	<input type="checkbox"/>	27-6430-02	Pre-designed 3 siRNA duplex standard purity, 2 nmol each	2 nmols
	<input type="checkbox"/>	27-6439-05	Pre-designed 3 siRNA duplex standard purity, 5 nmol each	5 nmols
	<input type="checkbox"/>	27-6430-10	Pre-designed 3 siRNA duplex standard purity, 10 nmol each	10 nmols



Research Use Only

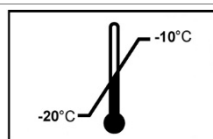

Storage

Store at -20°C to -10°C



Lot Number

Stated on product tube and packing slip



Expiry

Six Months from Date of Shipment



Instructions

Consult product manual



QR Code

Visit Gene Link website for product details

Handling & Storage

All siRNA are shipped lyophilized at room temperature. These are stable when kept lyophilized at room temperature. Follow established stringent RNase free handling conditions. The lyophilized siRNA duplex should be stored immediately at -20° C. The lyophilized siRNA is stable for ~6 months at -20° C.

*RNAi and siRNA

RNA interference (RNAi) is a specific and sequence dependent targeted gene silencing activity. RNAi acts by post transcriptional degradation of mRNA by small interfering RNAs (siRNA's) of the same sequence. The silencing approaches 100% and has to be empirically determined and optimized. Not every siRNA can effectively down regulate a gene. The process of RNA interference varies by individual siRNA while some do not exhibit any interference at all.

siRNA & RNAi Explorer™

RNAi Explorer™ from Gene Link is a series of product and services to aid researchers in exploring RNA interference. The online search and design algorithms for siRNA or shRNA are developed based on current known guidelines. You can place an order for the Guaranteed RNAi Explorer™ kit by simply providing Gene Link the appropriate Gene ID or Accession Number of your gene of interest. Alternatively, you can design your own siRNA or shRNA at Gene Link's website.

siRNA Synthesis

The siRNA duplexes provided are synthesized as a single strand oligo and annealed after complete deprotection and purification steps are completed. Gene Link utilizes the most reliable standard RNA synthesis chemistry. The RNA duplex is supplied lyophilized and ready to use.

Handling & Storage

Follow established stringent RNase free handling conditions. The lyophilized siRNA duplex should be stored immediately at -20° C. The lyophilized siRNA is stable for ~6 months at -20° C.

Guaranteed RNAi Explorer™ Kit

Please consult and save enclosed 'Order Specifications' sheet, this gives your gene specific siRNA information. This kit contains the following.

1. 3 siRNA duplexes 2 nmols each of 19+2 dT 3' overhangs. Supplied lyophilized, ready to dissolve and use. Please see order report for sequence information.
2. 1 negative control non-silencing siRNA duplex. 2 nmols supplied lyophilized, ready to dissolve and use.

Dissolve the lyophilized siRNA duplex in RNase-free deionized water as follows to give a final concentration of 20 μmolar or in any other desired concentration. Dissolve in an appropriate volume for any other quantity of siRNA.

siRNA Reconstitution				
	2 nmols	5 nmols	10 nmols	Work Sheet Space
RNase-free deionized water	100 μL	250 μL	500 μL	
Incubate the siRNA at 90° C for 1 minute, then at 37° C for 60 minutes. Aliquot in several tubes as required and store immediately at -20° C or preferably at -80° C.				

Guarantee: Gene Link guarantees a minimum of 70% silencing of your gene with at least one of the siRNA supplied. Gene Link will synthesize another set of 3 siRNA if evidence is provided that none of the provided siRNA was functional in achieving a minimum of 70% silencing.

Pre-designed 3 siRNA duplex standard purity

Please consult and save enclosed 'Order Specifications' sheet, this gives order specific data including the exact nmol quantity supplied of each duplex.

Custom siRNA Synthesis

Please consult and save enclosed 'Order Specifications' sheet, this gives order specific data including the exact nmol quantity supplied of each duplex.

Preparation of siRNA for Transfection

Dissolve the lyophilized siRNA duplex in RNase-free deionized water as follows to give a final concentration of 20 μ molar. Incubate the siRNA at 90° C for 1 minute, then at 37° C for 60 minutes.

Once it is dissolved, the siRNA should be aliquoted and stored at –80°C. For long term storage the siRNA can be ethanol precipitated in the following manner:

1. Add 1/10th volume of 10 M ammonium acetate.
2. Add 3 volumes of absolute ethanol and vortex briefly.
3. Store at –80°C.
4. When ready for use centrifuge at 14,000 x g for ten minutes.
5. Remove supernatant and leave the tube open to the air for 30 minutes in order to allow any residual alcohol to evaporate.
6. Dissolve in RNase-free deionized water as previously described.

Cell Transfection

1. Twenty-four hours prior to transfection, incubate a 24-well tissue culture plate with 2-10 x 10⁴ exponentially growing adherent cells per well.
2. Grow the cells overnight in an appropriate volume of growth medium. Cells should be approximately 40-70% confluent at the time of transfection.
3. Combine the siRNA and transfection reagent* according to the manufacturers specifications. In most cases this should be done in the absence of serum and antibiotics.
4. Remove the old medium from the cells and replace with fresh serum-containing medium.
5. Add the siRNA-transfection reagent complex.
6. Incubate the cells from 4-72 hours.
7. Perform appropriate assays to assess gene expression.

Partial list of commercially available transfection reagents: siPORT Amine Transfection Reagent from Ambion; BLOCK-IT Dicer RNAi Transfection kit from Invitrogen; TransIT-TKO siRNA Transfection Reagent from Mirus; RNAifect Transfection Reagent from Qiagen; GeneEraser siRNA Transfection Reagent from Stratagene and RiboJuice siRNA transfection Reagent from Novagen.

Controls for RNAi experiments

A negative control should also be included when performing siRNA experiments to eliminate the possibility of nonspecific silencing effects. For this purpose, Gene Link provides a siRNA sequence that has been shown to have no effect on gene silencing. The negative control siRNA is also available labeled with fluorescein, which allows transfection efficiency to be tracked by fluorescent microscopy. When monitoring transfection efficiency it should be noted that the pH indicator, phenol red, used in most culture media should be omitted as it interferes with fluorescent analysis.

Transfection conditions can be optimized further by the use of positive controls which consist of siRNA sequences that have been proven to be gene specific silencers.

Analyzing gene silencing

Gene silencing can be analyzed at either the protein or the mRNA level. Protein expression analysis is by western blotting, immunofluorescence, or FACS®. At the mRNA level gene silencing can be monitored by real-time RT-PCR, northern blotting, or microarray analysis. All gene expression data should be compared with levels of a non-related gene, such as beta actin or GAPDH, to normalize for variable amounts of RNA in different samples. Primers as well as Taqman probes and Molecular Beacons for beta actin and GAPDH are available from Gene Link.

Primer design for RT-PCR

Primers should be designed to amplify a region of the cDNA that spans an exon-exon boundary in order to prevent amplification from contaminating genomic DNA. Furthermore, primers should be designed to amplify the region flanking the siRNA binding site as partially degraded target mRNA, while no longer functional, may act as a template for PCR.

Troubleshooting & FAQ's

Low transfection efficiency

- ◆ When monitoring transfection efficiency by fluorescence uptake, measurements should be performed within 4-8 hours following transfection. After 8 hours the fluorescent signal may weaken even though the siRNA remains stable.
- ◆ Confirm that the medium used for formation of the siRNA-transfection reagent complex is serum-free.
- ◆ The transfection medium also should not contain polyanions such as heparin or dextran sulfate.
- ◆ Be sure the cell density at the time of transfection is optimal. The most favorable cell density for transfection is typically 40-70% confluence.
- ◆ Optimize the amount of transfection reagent by using a range of concentrations according to the manufacturer's recommendations.

Small or no gene silencing effect

- ◆ Monitoring of gene expression may be too soon to observe changes. This is particularly true if levels of protein expression are being monitored. A time course experiment should be performed to determine the optimal time for analysis.
- ◆ Perform a positive control experiment using a siRNA duplex that has been shown to be effective in gene silencing. Positive control siRNAs for human, mouse, and rat cell lines are available from Gene Link.

What are my siRNA delivery options?

Delivery of siRNA directly in cells can be achieved by using microinjection or electroporation. Another popular option is the use of transfection reagent. Specialized siRNA-delivery reagents are supplied by several companies. Please consult the transfection reagent vendor's protocol for detailed information for the exact requirements and procedure. Careful optimization of variable factors should be ensured for all initial transfection experiment. It is based on this and further optimization that reproducible gene knock out results will be obtained. Usually RNAi effect is seen within 4 hours and the maximum down regulation observed in 24-48 hrs. The effect lasts several cell generations and from 4-10 days depending on cell culture type.

How does purity of synthetic siRNA's affect RNAi?

RNAi is a sequence specific chain of events. Chemical synthesis of siRNA's is based on coupling of bases to yield a particular sequence. The yield and purity depends on the coupling efficiency. Gene Link siRNA's can be used without further purification, but Gene Link recommends purified siRNA's for use in transfection.

What concentration of siRNA is most effective?

As low as 1 nM concentration of siRNA have been shown to be effective in exhibiting RNAi. Initial experiments should be done at varying concentrations from 1-10 nM. Some reports have used as high as 25nM concentrations. High quality siRNA's should be used.

How do I quantify down regulation?

RNAi down regulates a gene function without actually interacting with the gene. The subtle action is by mRNA degradation. Thus the degree of RNA interference achieved is directly proportional to the level of mature mRNA and the translated proteins. The options are:

1. Measurements of target protein (enzyme) activity. This option is suitable if a robust assay is available or has been in prior use. The assay would vary by the nature of the protein product.
2. Measurement of target mRNA level. This is the preferred method as it directly quantifies the level of mRNA. Quantitative PCR is very effective in measuring relative amount of target sequence. This can be achieved simply by SYBR green or by the use of TaqMan or Molecular Beacons.

References for further information

1. Elbashir, S.M., Harborth, J., Lendeckel, W., Yalcin, A., Weber, K. and Tuschl, T. (2001) Nature, 411, 494-498 .
2. Yu Jenn-Yah, DeRuiter, S.L, and Turner, D.L. (2002) PNAS, 99, 6047-6052.
3. Gunther, S., Hanna, D., William B., K., Nicole, R., Marc H., S., Clare M., W., Sandra L., S. and M.A., S. (2004) Mol. Biol of the Cell , 15, 256-267.
4. Yang H., H., Soung H., J. and Seung H., Y. (2003) The Journal of Biological Chemistry, 278, 40581-40589
5. Jens, H., Sayda M., E., Kim, B., Thomas, T. and Klaus, W. (2001) Journal of Cell Science, 114, 4557-4565
6. Hannon, G.J. (2002) RNA interference. Nature, 418, 244-251
- 7.. Fire, A., Xu, S., Montgomery, M.K., Kostas, S.A., Driver, S.E. and Mello, C.C. (1998) Nature, 391, 806-811.



***RNAi and siRNA**

RNA interference (RNAi) is a specific and sequence dependent targeted gene silencing activity. RNAi acts by post transcriptional degradation of mRNA by small interfering RNAs (siRNA's) of the same sequence. The silencing approaches 100% and has to be empirically determined and optimized. Not every siRNA can effectively down regulate a gene. The process of RNA interference varies by individual siRNA while some do not exhibit any interference at all.



Related Products Ordering Information

DNA & RNA Precipitation Solutions

Product	Catalog No.	Unit Size
DNA & RNA Precipitation Solutions Pack (contains the following; Glycogen Solution 10 mg/mL; 1 mL [40-5112-01]; Linear Acrylamide Solution 5mg/mL; 1 mL [40-5113-01] LiCl RNA Precipitation Solution [40-5131-05]; Sodium Acetate DNA & RNA Precipitation Solution [40-5132-05]; Sodium Chloride DNA & RNA Precipitation [40-5134-05] and Ammonium Acetate 7.5M DNA & RNA Precipitation Solution [40-4135-05])	40-5130-00	1 Pack
Glycogen Solution 10 mg/mL; 1 mL	40-5112-01	1 mL
Linear Acrylamide Solution (Linear polyacrylamide, LPA; 5mg/mL); 1 mL	40-5113-01	1 mL
LiCl RNA Precipitation Solution (7.5M LiCl, 50 mM EDTA pH 8.0); 50 mL	40-5131-05	50 mL
Sodium Acetate 3M pH 5.5; DNA & RNA Precipitation Solution; 50 mL	40-5132-05	50 mL
Potassium Acetate 3M pH 5.5; DNA & RNA Precipitation Solution; 50 mL	40-5133-05	50 mL
Sodium Chloride 5M DNA & RNA Precipitation; 50 mL	40-5134-05	50 mL
Ammonium Acetate 7.5M DNA & RNA Precipitation Solution; 50 mL	40-5135-05	50 mL
Ammonium Acetate 5 M DNA & RNA Precipitation Solution; 50 mL	40-5136-05	50 mL

Related Products Ordering Information

DNA & RNA Reconstitution Solutions

Product	Catalog No.	Unit Size
DNA & RNA Reconstitution Solutions Pack (contains 50 mL each of DEPC Treated Water [40-3000-05], Nuclease Free Water (DEPC Free) [40-3001-05], TE pH 7.0 [40-5011-05] and RNA Reconstitution Solution [40-5014-05])	40-3000-00	1 Pack
RNA Reconstitution & Storage Solution (1 mM Sodium Citrate pH 6.4); 10 X 1.6 mL	40-5014-16	10 X 1.6 mL
RNA Reconstitution & Storage Solution (1 mM Sodium Citrate pH 6.4); 50 mL	40-5014-05	50 mL
TE Buffer 1X solution pH 7.0; 50 mL	40-5011-05	50 mL
TE Buffer 1X solution pH 7.5; 50 mL	40-5012-05	50 mL
TE Buffer 1X solution pH 8.0; 50 mL	40-5013-05	50 mL
Nuclease Free Water (DEPC Free); 10 X 1.6 mL	40-3001-16	10 X 1.6 mL
Nuclease Free Water (DEPC Free); 50 mL	40-3001-05	50 mL
Nuclease Free Water (DEPC Free); 500 mL	40-3001-50	500 mL
Nuclease Free Water (DEPC Free); 1L	40-3001-01	1 L
DEPC Treated Water; 10 X 1.6 mL	40-3000-16	10 X 1.6 mL
DEPC Treated Water; 50 mL	40-3000-05	50 mL
DEPC Treated Water; 500 mL	40-3000-50	500 mL
DEPC Treated Water; 1L	40-3000-01	1 L

Document Warranty and Liability

Information in this document is subject to change without notice. This document and all information presented in this document are written as a guide. Gene Link, Inc. does not warrant this document to be free of errors and assumes no responsibility for any errors that may appear in this document.

Gene Link disclaims all warranties with respect to this document, expressed or implied, including but not limited to those of merchantability or fitness for a particular purpose. In no event shall Gene Link be liable, whether in contract, tort, warranty, or under any statute or on any other basis for special, incidental, indirect, punitive, multiple or consequential damages in connection with or arising from this document, including but not limited to the use thereof.

Website

As the receipt of information on the Internet is highly dependent upon factors, including without limitations to, the user's computer, browser, operation system, etc., information may be perceived incorrectly. Therefore, Gene Link does not warrant or guarantee that the information contained on its website www.genelink.com is error free.

Product Warranty and Liability

Warranty: Gene Link makes no warranty of any kind, specifically disclaims and excludes all other warranties of any kind or nature, directly or indirectly, express or implied, including, without limitation, as to the suitability, productivity, durability, fitness for a particular purpose or use, merchantability, condition, or any other matter with respect to Gene Link products. Gene Link products are for research purposes only including custom products. There is no warranty or claim of its performance for any specific research application. All Gene Link products are guaranteed to meet or exceed the specifications stated. Each Gene Link product is shipped with documentation stating specifications and other technical information. If the product fails to meet the stated specifications the sole remedy is prompt replacement by Gene Link or within 30 days of purchase a refund of the purchased price.

Liability: Under no circumstances shall Gene Link be liable for any damages directly or indirectly related to Gene Link's products and services. Whether direct, incidental, foreseeable, consequential, or special (including but not limited to loss of use, revenue or profit), whether based upon warranty, contract, tort (including negligence) or strict liability arising in connection with the sale or the failure of Gene Link products to perform in accordance with the stated specifications.

Research Use Only. Not for use in diagnostic or clinical procedures.

Notice to Purchaser: The purchase of this product conveys to the purchaser the limited, non-transferable right to use the purchased amount of the product only to perform internal research for the sole benefit of the purchaser. No right to resell this product or any of its components is conveyed expressly, by implication, or by estoppel. This product is for internal research purposes only and is not for use in commercial applications of any kind, including, without limitation, quality control and commercial services such as reporting the results of purchaser's activities for a fee or other form of consideration. For information on obtaining additional rights, please contact support@genelink.com.

© 2015 Gene Link Inc. All rights reserved.

The trademarks mentioned herein are the property of their respective owners.

Gene Link, Inc.
190 Saw Mill River Road
Hawthorne, NY 10532
USA

Tel: (914) 769-1192
Email: support@genelink.com
www.genelink.com