



Product Specifications


Custom Oligo Synthesis, antisense oligos, RNA oligos, chimeric oligos, Fluorescent dyes, Affinity Ligands, Spacers & Linkers, Duplex Stabilizers, Minor bases, labeled oligos, Molecular Beacons, siRNA, phosphonates Locked Nucleic Acids (LNA); 2'-5' linked Oligos

Oligo Modifications

For research use only. Not for use in diagnostic procedures for clinical purposes.

DBCO-dT

| | |
|--------------------------|-----------------|
| Category | Click Chemistry |
| Modification Code | DBCO-dT |
| Reference Catalog Number | 26-6927 |
| 5 Prime | Y |
| 3 Prime | N |
| Internal | N |
| Molecular Weight(mw) | 773.77 |

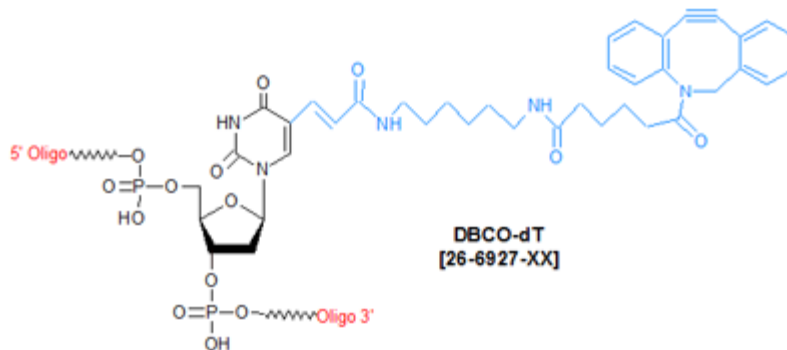

[Click here for a complete list of Click Chemistry Oligo Modifications](#)

DBCO-TEG, DBCO-dT and DBCO Serinol are discontinued modifications due to their instability during direct oligo synthesis protocols. We offer a range of post synthesis DBCO NHS modifications as alternate. See above related modifications.

DBCO (dibenzocyclooctynes, DBCO) conjugation chemistry is based on the reaction of a dibenzylcyclooctyne (DBCO) linker with an azide linker to form a stable triazole. The dibenzocyclooctyne group (DBCO) allows Copper-free Click Chemistry to be done with live cells, whole organisms, and non-living samples. DBCO groups will preferentially and spontaneously label molecules containing azide groups (N₃). Within physiological temperature and pH ranges, the DBCO group does not react with amines or hydroxyls, which are naturally present in many biomolecules. Reaction of the DBCO group with the azide group is significantly faster than with the sulfhydryl group (SH, thiol).

Cyclooctyne-based modifications offers the ease of copper-free click reagents. These are simple to use and has excellent click performance in 17 hours or less at room temperature. Gene Link offers 5' DBCO-TEG for preparing oligos with 5' DBCO and a 15 tom triethylene glycol spacer arm, **DBCO-dT is used for inserting a DBCO group preferably at the 5' or internally within 3-4 bases from the 5'** and DBCO-sulfo-NHS Ester is also offered for post-synthesis conjugation reactions. DBCO-modified oligos may be conjugated with azides in organic solvents, such as DMSO, or aqueous buffers. Depending on the azide used, the reaction will go to completion in 4-17 hours at room temperature.

Glen Report 27.1: Technical Brief - DBCO-dT - An Unusual Case of Iodine Sensitivity